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


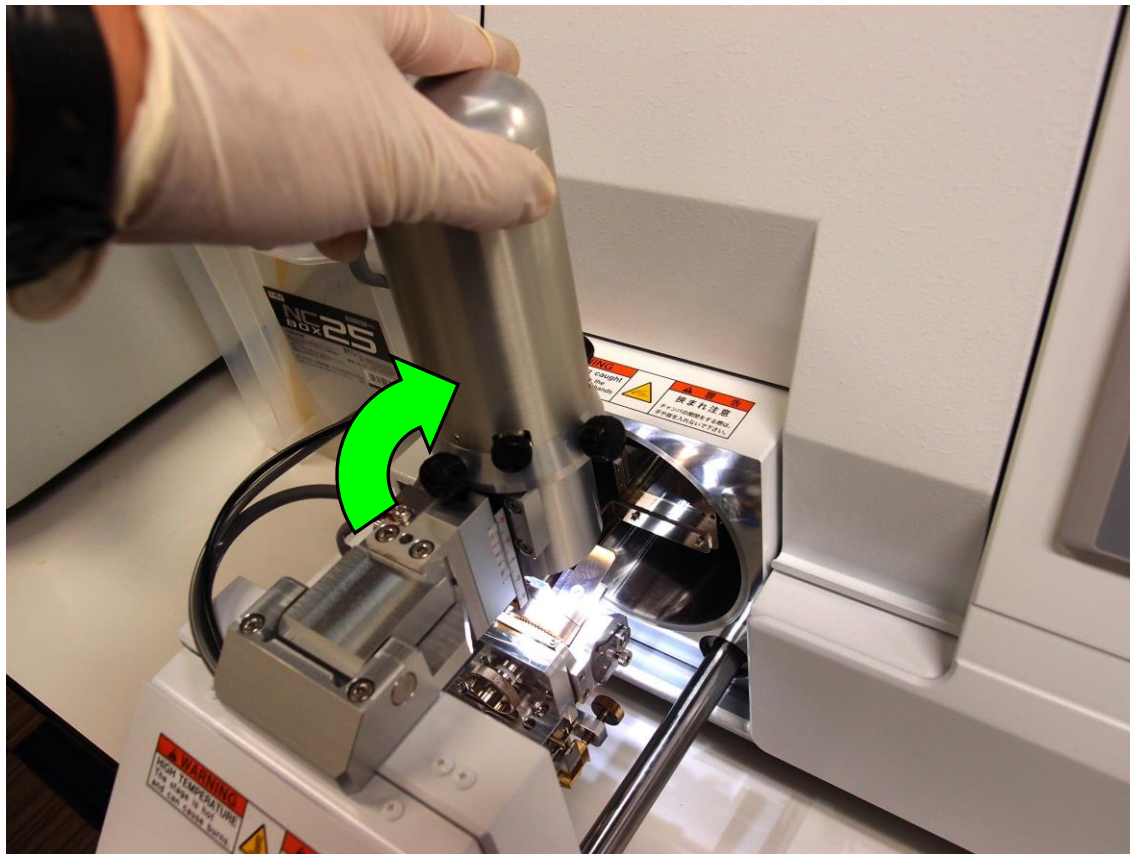
The image shows a control panel interface with a blue background. At the top, there are five menu buttons: 操作画面, プリセット, ガス最適化, カメラ, and 設定. A green arrow points to the 'カメラ' button. Below the menu buttons, there are several control elements:

- On the left, there are two circular buttons labeled 'VENT' and 'EVAC', and a vertical black bar with 'Pa' below it.
- In the center, there are three stacked boxes: the top one shows '5.0 kV', the middle one shows 'Arガス流量 3.3', and the bottom one shows 'イオンビーム On OFF'.
- To the right of these boxes, there is a control panel with a '0  $\mu A$ ' display and a 'タイマー 04:00' display. Below these is a large '00:00' display.
- On the far right, there are four stacked buttons: '自動開始モード', '自動電源停止モード', '試料回転ホルダー', and '仕上げ加工モード'. At the bottom right is a large button with a play icon and the text '開始'.

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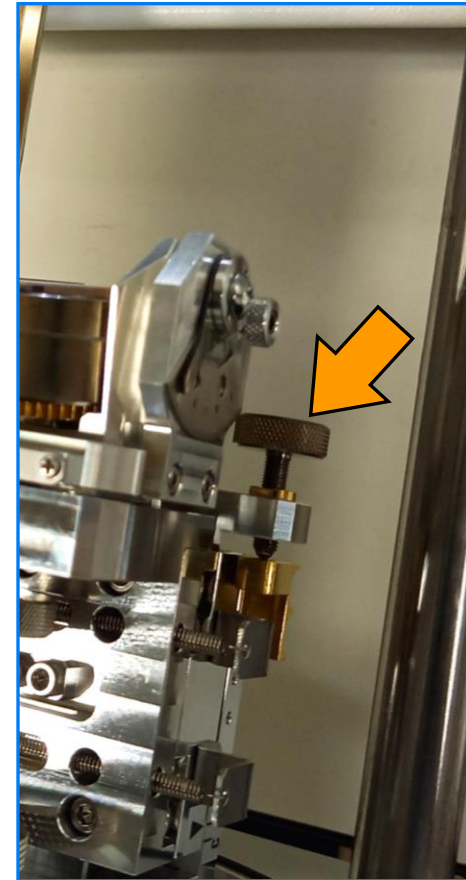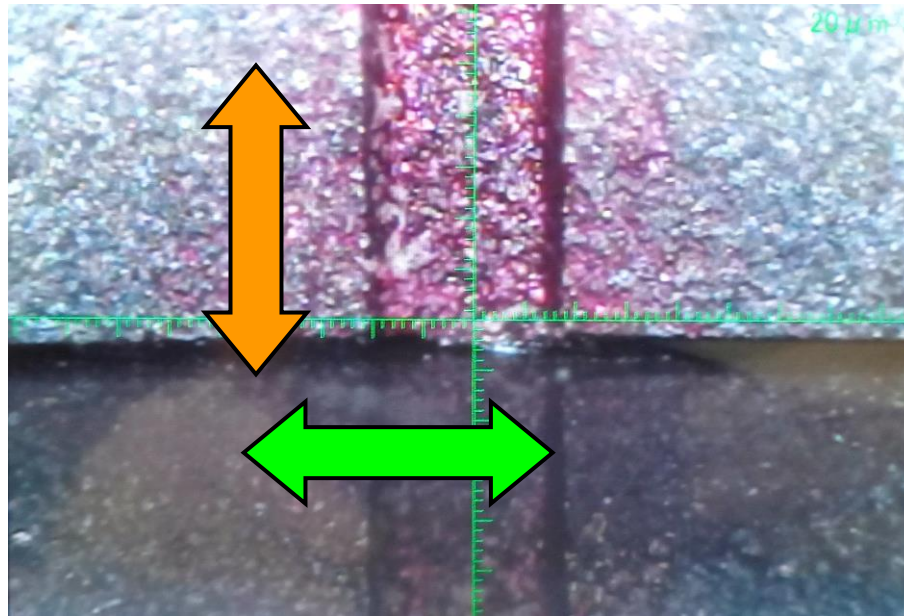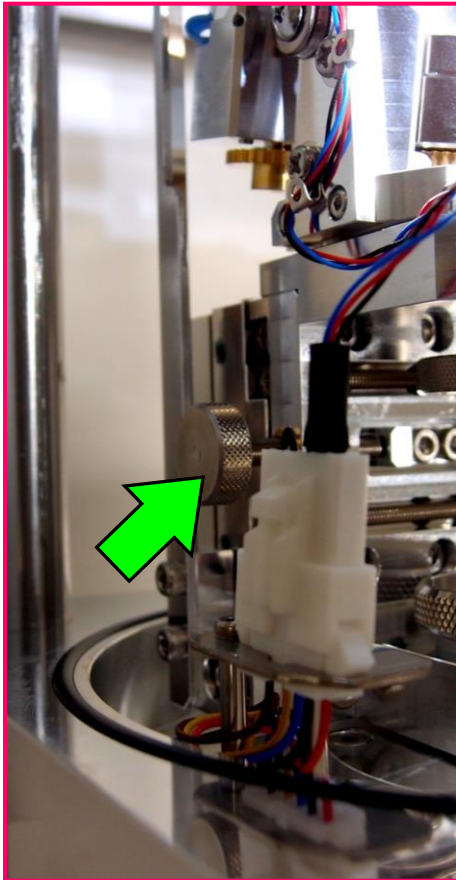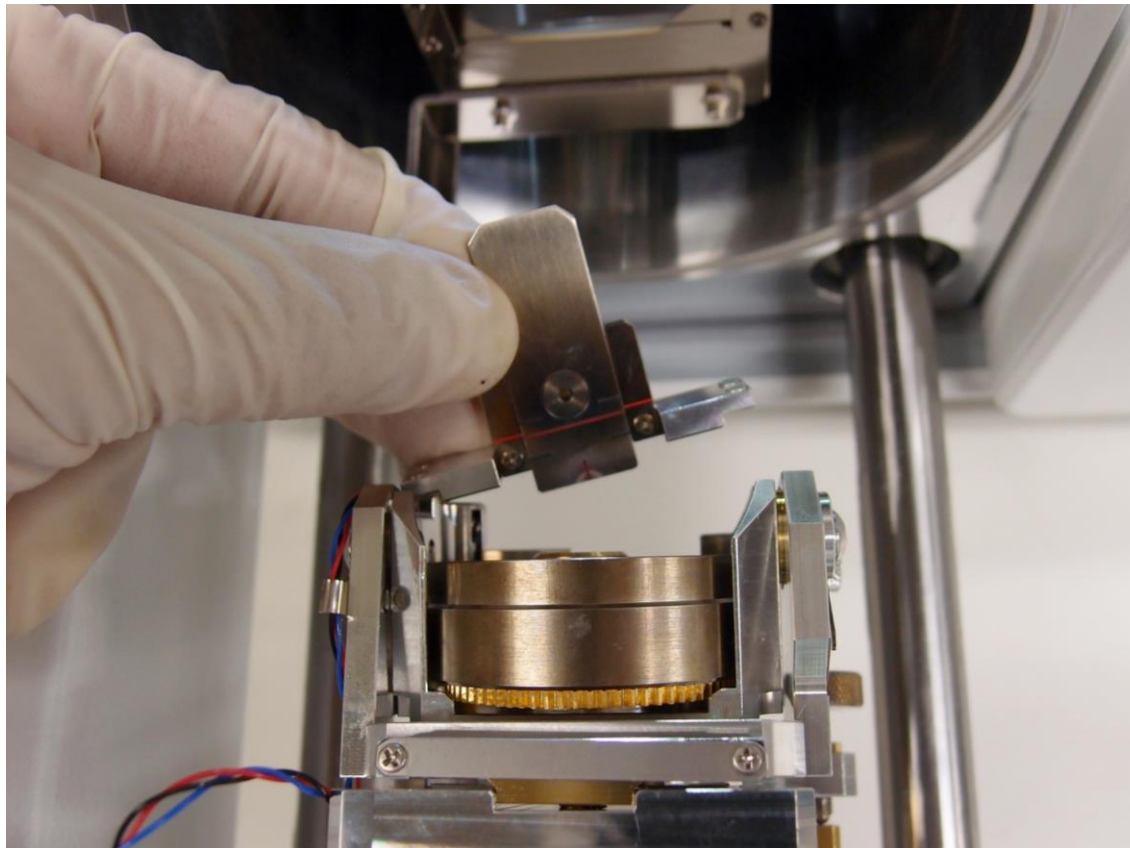
The control panel features a top navigation bar with five tabs: 操作画面, プリセット, ガス最適化, カメラ, and 設定. On the left side, there are two circular buttons: VENT (white) and EVAC (black with white text). A green arrow points to the EVAC button. Below these buttons is a vertical black bar with the label Pa at the bottom left. The main display area contains several data fields: a large display showing 5.0 kV, a display showing 0 μA with up and down arrow controls, a display showing Arガス流量 3.3, a display showing イオンビーム On OFF, and a large digital timer showing 00:00 with a label タイマー 04:00 above it. On the right side, there are five rectangular buttons: 自動開始モード, 自動電源停止モード, 試料回転ホルダー (highlighted in black with green text), 仕上げ加工モード, and 開始 (with a play button icon).

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